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Table 1-3. Operating Specifications for Model 05-STP Series (cont'd)

Parameter	Minimum	Typical	Maximum	Units
Shock				
Impulse (IEC68.2.27)		30x11		gxmsec
Vibration (IEC68.2.6)	10		55	Hz
Warm-up		15	40	minutes

Optical				
Model 05-STP-910 Single Frequency Laser System				
Parameter	Minimum	Typical	Maximum	Units
Output Range				
Power (633 nm), Typical	0.35		0.70	mW
Frequency ($\nu - \nu_{o(1)}$)	400		600	MHz
Beam Diameter ($1/e^2$)		0.48		mm
Beam Divergence (Full Angle)		1.70		mrاد
Spatial Mode (TEM_{00})	100			%
Optical Noise			0.10	%RMS
Polarization	1000:1	5000:1		
Power Stability				
1 Hour		± 0.1		%
8 Hour		± 0.1		%
1 month		± 0.1	± 0.2	%
Tempco			0.18	%/ $^{\circ}C$
Frequency Stability ¹				
1 Hour		$\pm$		MHz
8 Hour		$\pm$		MHz
1 Month			± 0	MHz
Tempco		3.60		MHz/ $^{\circ}C$

Notes:

1. See Section 1.7.

Table 1-3. Operating Specifications for Model 05-STP Series (cont'd)

Optical				
Model 05-STP-912 Single Frequency Laser System				
Parameter	Minimum	Typical	Maximum	Units
Output Range				
Power (633 nm), Typical	0.60		1.40	mW
Frequency ($\nu - \nu_{o(1)}$)	50		600	MHz
Beam Diameter ($1/e^2$)		0.54		mm
Beam Divergence (Full Angle)		1.50		mrاد
Spatial Mode (TEM_{00})	100			%
Optical Noise			0.10	%RMS
Polarization	1000:1	5000:1		
Power Stability				
1 Hour		±0.1		%
8 Hour		±0.1		%
1 month		±0.1	±0.2	%
Tempco			0.18	%/°C
Frequency Stability ¹				
1Hour		±		MHz
8 Hour		±		MHz
1 Month			±0	MHz
Tempco		3.60		MHz/°C

Notes:

1. See Section 1.7.

05-STP-910/912

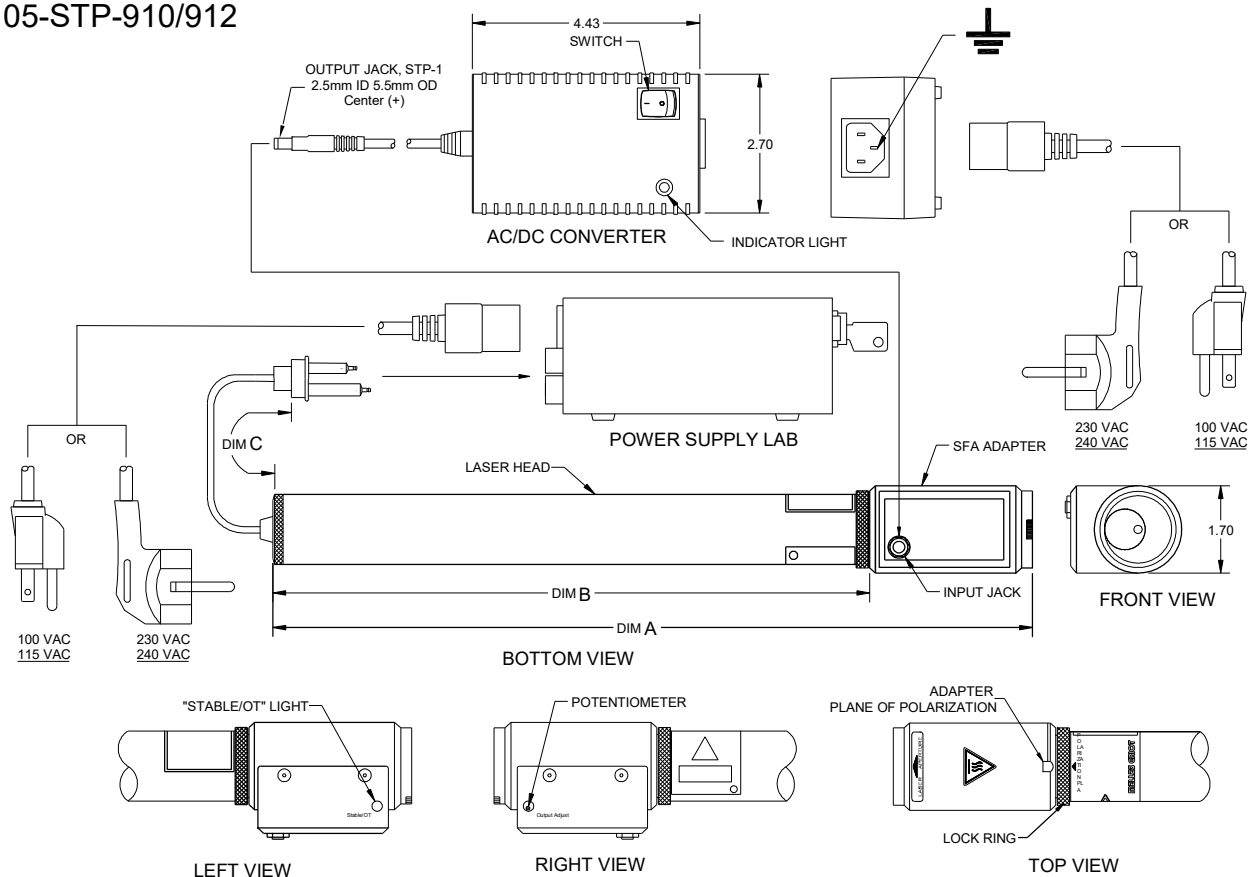
Figure 1-4

MODEL	DIM A	DIM B	DIM C
05-STP-910-XXX	10.20 (259.1)	6.96 (176.8)	72.00 (1828.8)
05-STP-912-XXX	11.89 (302.0)	8.65 (219.7)	72.00 (1828.8)

Figure 1-4. Mechanical Specifications of Models 05-STP-910 and 05-STP-912

1.7. Doppler Output Power/Frequency Function

Understanding the basic properties of a specific laser allows the output frequency of the Models 05-STP-910 or 05-STP-912 system to be accurately tuned over a large portion of the available doppler gain profile.

The output power of the Model 05-STP-910 or 05-STP-912 system as a function of output frequency (ν) can be described by the following relationship.

Formula 1

$$P_o(\nu) = A(n^2 \exp[-B(\nu - \nu_o)^2 / \nu_o^2])$$

where

- * P_o is the output power in mW
- * ν is the absolute output frequency
- * n is the tube excitation parameter (total single pass gain/total losses)
- * A, B are coefficients that describe various physical constants and tube design parameters
- * ν_o is the transition frequency center-line that is established as 473.6122×10^{12} Hz.

The "Doppler" output profile can be described in terms of output frequency with reference to the absolute transition frequency due to the magnitude of values encountered in the visible spectrum. In this case formula 1 is simply rewritten as:

$$P_o(\lambda) = A(n^2 \exp[-B\lambda^2 / \nu_o^2])$$

Figure 1-7 is a theoretical plot of the above equation versus λ for two conditions of n . Curve 1 depicts a normal condition of n that exists for a tube at the beginning of its operating life.

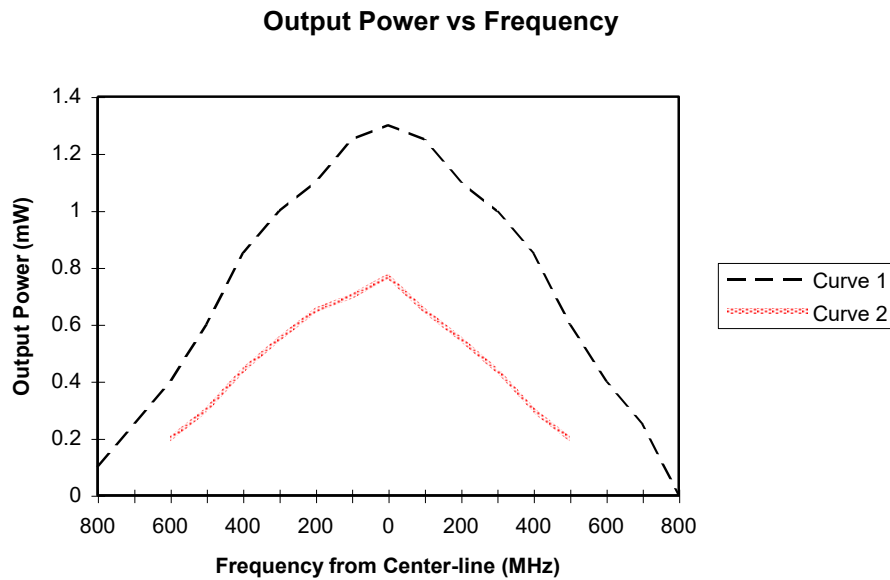
Figure 1-5. Output Power vs. Frequency

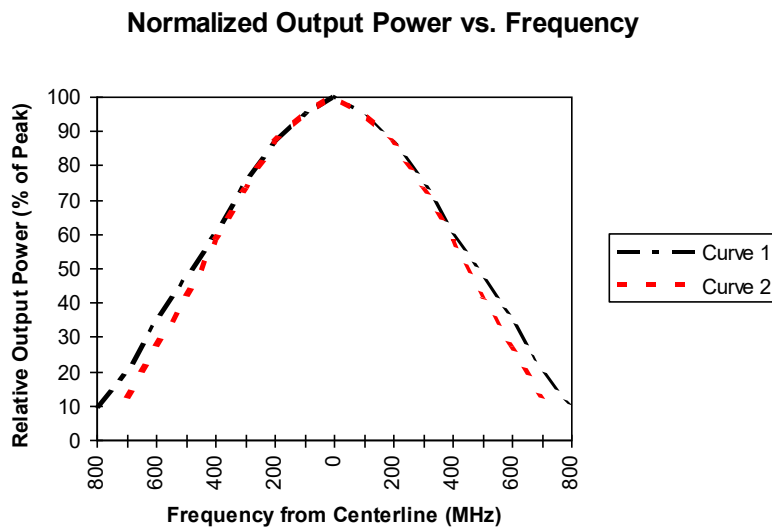
Figure 1-6. Normalized Output Power vs. Frequency

Curve 2 depicts the doppler profile that exists at the end of the operating life if n falls to 80% of its original value. This gain decrease may be from gas cleanup or increased absorption loss changes of the cavity mirrors.

The peak output power available (at center-line) drops to nearly half its original value.

If one normalizes both curves in Figure 1-7, it can be noted that this extreme degradation does not appreciably effect the shape of the doppler curve. Thus, if one were to know the peak value of output power (at center-line), a reasonably accurate calibration of output frequency could be attained by using a power meter a calibration curve similar to Figure 1-9.

1.8. Frequency Calibration

As explained in Section 1.7, the Model 05-STP-910, 05-STP-912, single frequency laser system may be tuned over a wide range of frequencies from center-line. Using a calibration curve, such as the one shown in Figure 1-9, the output frequency can be calibrated to typically better than $\pm 0\%$ accuracy of a given frequency on the "Blue" side of the doppler output profile for the 632.991 nm transition in Ne₂₀.

Figure 1-9 was constructed from measurements of a typical model 05-STP-910 system producing peak center-line power of about 1.3mW. As shown in Figure 1-8, the shape of this curve will not deviate substantially over the useful life of the laser.

A laser radiometer may be used for calibration by noting the peak center-line power during the warm-up period while in the "Peak" detection mode. The desired frequency can then be adjusted by looking up on the ordinate of Figure 1-9, the appropriate percentage of peak center-line power for the desired frequency. By multiplying this percentage by the peak level noted on the power meter, the output power may be adjusted for the desired frequency.

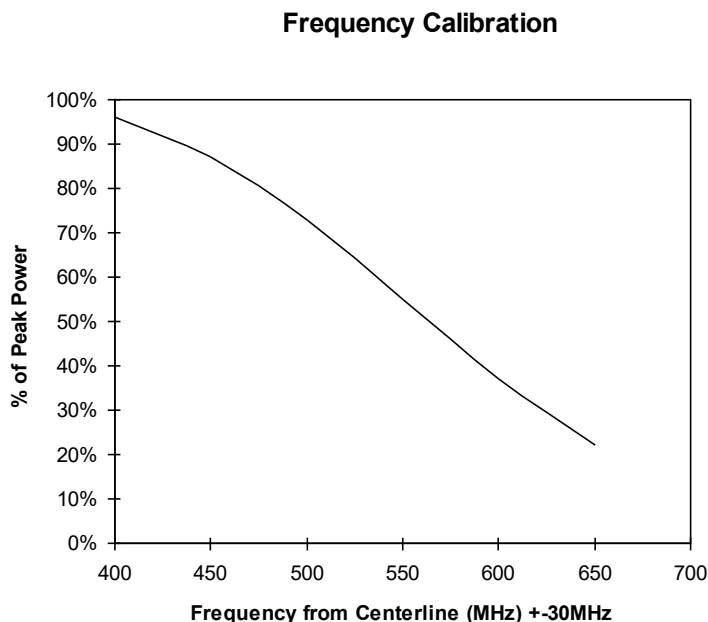
Figure 1-7. Calibration Curve

APPENDIX A: WARRANTY

Pacific Lasertec (“Manufacturer”) warrants its products to be free from defects in materials or workmanship for a period of one year from the date of shipment from the factory. The Manufacturer’s liability is limited to replacing, repairing or issuing credit, in our sole discretion, for any products that are returned by the original purchaser during the warranty period and fail to meet specifications. The Manufacturer makes no warranty that the products are fit for the use or purpose to which they may be put by the buyer, whether or not such use or purpose has been disclosed to the Manufacturer in specifications or drawings previously or subsequently provided, or whether or not the products are specifically designed and/or manufactured for buyer’s use or purpose. The Manufacturer’s liability for any claim for loss or damage arising out of the sale, resale or use of any of its products shall in no event exceed the selling price of the unit. Please contact the Manufacturer for any further required information:

Pacific Lasertec
215 Bingham Drive, Suite 110
San Marcos, California 92069, U.S.A.
Phone: (760) 539-7169

Email: contact@pacifclasertec.com

Please visit: <https://pacifclasertec.com/contact>
for International Distributors and Representatives

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© 2019 Pacific Lasertec